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The screenshot displays the EGCloud registration interface. At the top, the EGCloud logo is centered. Below it, a progress bar indicates three steps: 1. 账户信息 (Account Information), 2. 辅助信息 (Auxiliary Information), and 3. 注册完成 (Registration Complete). The current step, '辅助信息', contains several input fields: '姓名' (Name) with the placeholder '姓名', '性别' (Gender) with radio buttons for '男' (Male) and '女' (Female), '公司邮箱' (Company Email) with the placeholder 'info@company.com', '主页' (Homepage) with the placeholder 'https://www.google.com', and '备注' (Remarks) with the placeholder '请输入备注'. A button labeled '注册' (Register) and a button labeled '上一步' (Previous Step) are positioned below the fields. A link for '使用已有账户登录' (Log in with existing account) is located below the buttons. A grey button labeled '省略部分项目...' (Omit some items...) is also visible. At the bottom of the page, the copyright notice 'Copyright © 2019 EGCloud 出品' is displayed.

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The screenshot shows the EGCloud web interface for managing experimental sites. The top navigation bar is blue and contains the logo, a menu icon, the text '欢迎进入 EGCloud', a user profile icon, '欢迎您, 系统用户', and a '退出登录' button. The left sidebar has a menu with '首页', '实验场所', '设备登录', '实验管理', '代理设置', and '个人中心'. The '实验场所' menu item is selected and highlighted in blue. The main content area is titled '实验场所' and features a search bar with the placeholder '请输入实验场所名称', a '查询' button, and a '重置' button. Below the search bar is a '+ 新增' button. A selection bar indicates '已选择 0项' and a '清空' button. A table displays the following data:

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At the bottom right of the table area, there is a pagination control showing '1-2 共2条' and a dropdown menu for '10条/页'. The footer of the page contains the text 'Copyright © 2019 EGCloud'.

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The screenshot displays the '个人中心' (Personal Center) page in the EGCloud system. The page is divided into a left sidebar and a main content area. The sidebar contains navigation links: 首页 (Home), 实验场所 (Experiment Location), 设备登录 (Device Login), 代理设置 (Proxy Settings), 实验编辑 (Experiment Edit), 实验一览 (Experiment Overview), 历史记录 (History), and 个人中心 (Personal Center). The main content area is titled '系统用户' (System User) and features a profile picture and a '标签' (Tags) section with 'system expert' and '+New Tag'. The '基本设置' (Basic Settings) tab is selected, showing fields for '姓名' (Name) with value '系统用户', '性别' (Gender) with radio buttons for '男' (Male) and '女' (Female), '联系电话' (Contact Number) with value '185666662', '简介' (Introduction) with a text area, and '电子邮件' (Email) with value 'username@example.com'. A '保存' (Save) button is located at the bottom of the form. The footer of the page reads 'Copyright © 2019 EGCloud'.

END

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Copyright © 2019 EGCLOUD